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Installation Manual for RS232/RS422-RS485 Interface Converter

Content

6.	Service	5
1.	Information notice regarding the manual	6
2.	Intended use	6
3.	Unit description	7
4.	Scope of delivery of RS232/RS422-RS485 interface converter kit	8
5.	Networking	8
5.1.	Bus termination / RS485 operation	9
6.	Service	9

1. Information notice regarding the manual

This installation manual contains a short description how to connect and install the RS232/RS422-RS485 interface converter and the drivers needed. It is expected that the Interface Converter has been purchased at the BINDER GmbH and thus the correct internal settings have already been carried out. It is also assumed that only devices and cables of the BINDER GmbH offered specially for this purpose will be used. To allow also different kinds of use offered by the RS232/RS422-RS485 interface converter, we include next to this manual also the original operation manual of the producer W & T.

2. Intended use

The RS232/RS422-RS485 interface converter is suitable to network temperature chambers with RS422 or RS485 interface, measurement devices, and the Alarm Box AB01 of the BINDER GmbH via the RS232 computer interface.

At any computer the communication software APT-COM™ of the BINDER GmbH or an existing data acquisition system can be used to control the temperature chambers connected to the RS422 network or RS485 network and to receive and save the chamber measurement data.

3. Unit description

The RS232/RS422-RS485 interface converter is independent electrical equipment with external power supply via a 24V AC power supply unit. It can be loosely installed or mounted to clamping profile bars in distributor boxes. By internal DIP switches and plug bridges different operational modes can be set. The settings necessary for operating temperature chambers of the BINDER GmbH have already been pre-set by the BINDER GmbH.

The operating mode for BINDER RS422 chambers according to the original W&T user manual is "RS422, RS 485, 4-wire-bus-master". The operating mode for BINDER RS485 chambers according to the original W&T user manual is "RS485, 2-wire without echo, with automatic control". For standard settings for BINDER chambers refer to Figure 3.

The connection to a RS232 interface is established using a Zero-Modem cable. The connection to a temperature chamber of the BINDER GmbH or to a plug distributor for connecting several temperature chambers is established using a 9-poles Sub-D socket and a special cable. Both connection cables and the power supply unit are included with the scope of delivery.

The connected temperature chamber(s) are addressed via this COM interface and the address that can be individually adjusted at each chamber controller.

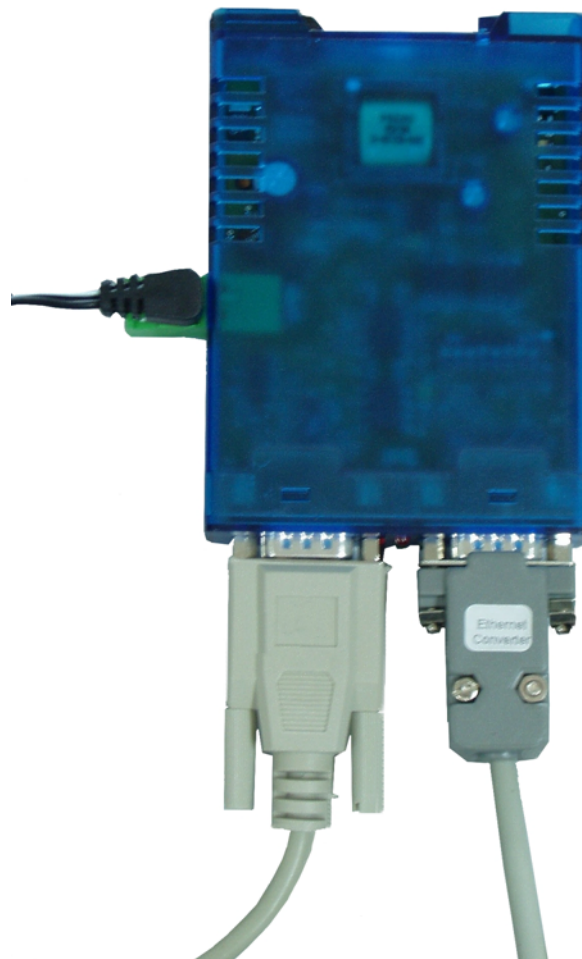
Figure 1: Connected interface converter

4. Scope of delivery of RS232/RS422-RS485 interface converter kit

Complete kit

Description	Art.no.
RS232/RS422-RS485 interface converter kit	8012-0556

Components:

Description	Art.no.
RS232/RS422-RS485 interface converter with original description of producer W & T (only needed in case of problems during start-up)	5021-0022
Zero-Modem cable 3 m for connection to RS232 interface	5023-0115
Connection cable 5m for connecting the interface converter to an individual chamber or to a plug distributor RS422 from BINDER	5023-0095
Manual for connecting the RS232/RS422-RS485 interface converter	7001-0101

Option:

Description	Art.no.
Extension cable 50m for connecting the connection cable 5m to an individual chamber or to a plug distributor RS422 from BINDER	5023-0117

5. Networking

All temperature chambers are connected via the same RS232/RS422 interface converter.

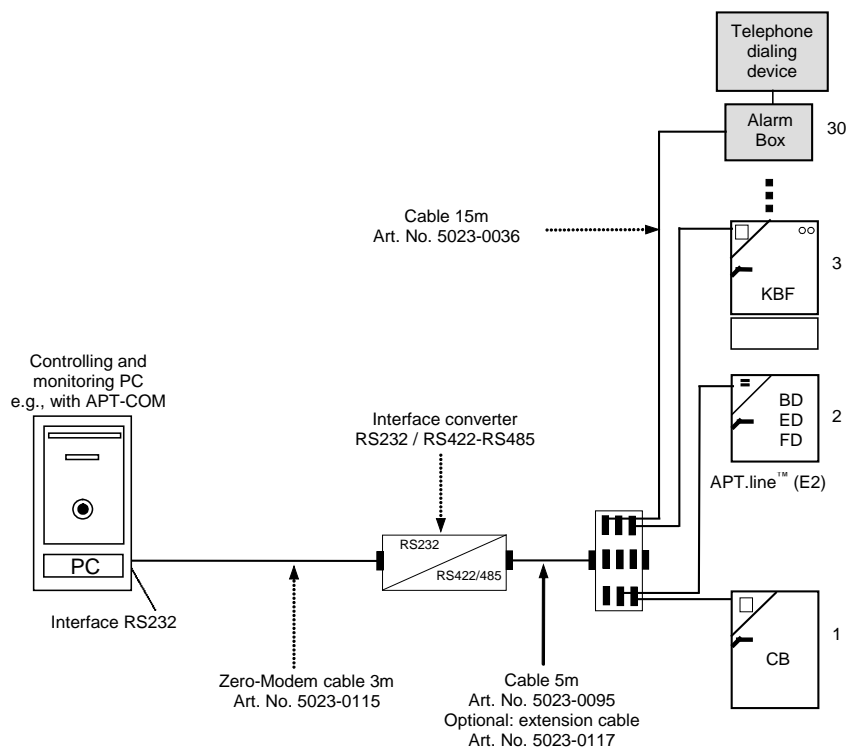
Figure 2: Networking via RS232/RS422-RS485 interface converter

The optional extension cable 50m (5023-0117) can be connected behind the 5m cable (5023-0095).

The RS232/RS422-RS485 interface converter is now completely installed. The temperature chambers are now registered in the communication software APT-COM™ with the COM number that has been assigned to the RS232/RS422-RS485 interface converter they are altogether connected to. The chamber address required for the unambiguous chamber identification must be entered by hand individually and only once at the controller of each temperature chamber or measurement device. For a description see the respective operating manual of the chamber or device bearing an RS422 interface or RS485 interface.

! To integrate RS485 chambers into an existing RS422 network, please use the RS485/RS422 interface converter:

230V version: Art. no. 8012-0589, 115V version: Art. no. 8012-0599

5.1. Bus termination / RS485 operation

For the connection of RS485 networks, DIP switches 1 and 2 must be set to ON

To connect temperature chambers from BINDER, the recommended operating mode of the interface converter is termination enabled (BINDER factory setting).

Some controllers, like RD3, MB1, MB2, have an internal termination. When combining different controllers, communication errors can occur which are indicated by APT-COM™ as „timeout” or “CRC error”. In this case it can be necessary to disable termination at the interface converter.

To disable termination, set DIP switches 6 and 7 in the interface converter to Off. You can safely try out the correct setting. No damage of controller or interface converter will occur.

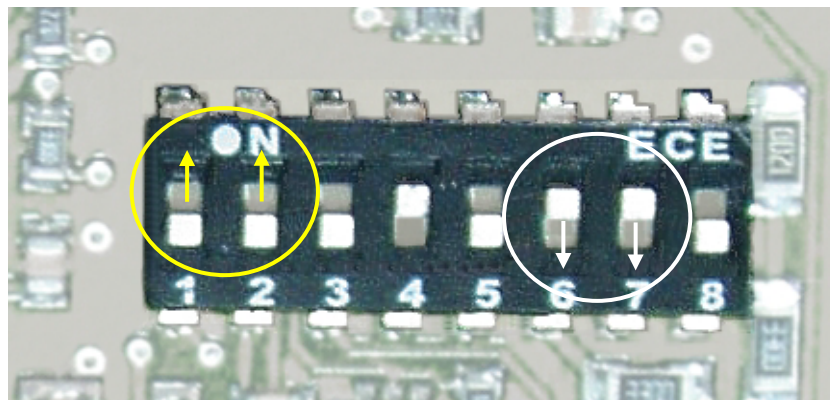
Figure 3: Termination setting via DIP switches 6 and 7

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