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Installation Manual for RS485/RS422 Interface Converter

Content

1.	Information notice regarding the manual	7
2.	Intended use	7
3.	Unit description	8
4.	Scope of delivery of RS485/RS422 interface converter kit	9
5.	Networking	10
5.1.	Bus termination	11
6.	Service	11

1. Information notice regarding the manual

This installation manual contains a short description how to connect and install the RS485/RS422 interface converter. It is expected that the Interface Converter has been purchased at the BINDER GmbH and thus the correct internal settings have already been carried out. It is also assumed that only devices and cables of the BINDER GmbH offered specially for this purpose will be used. To allow also different kinds of use offered by the RS485/RS422 interface converter, we include next to this manual also the original operation manual of the producer W & T.

2. Intended use

The RS485/RS422 interface converter is suitable to integrate temperature chambers with a RS485 interface into an existing RS422 network.

3. Unit description

The RS485/RS422 interface converter is independent electrical equipment with external power supply via a 24V AC power supply unit. It can be loosely installed or mounted to clamping profile bars in distributor boxes. By internal DIP switches and plug bridges different operational modes can be set. The settings necessary for operating temperature chambers of the BINDER GmbH have already been pre-set by the BINDER GmbH.

The operating mode for BINDER chambers according to the original W&T user manual is on the RS422 connection side "RS 485, 4-wire-bus-master" and on the RS485 connection side "RS485, 2-wire-bus-master". For standard settings for BINDER chambers refer to Figure 3.

Both connection cables and the power supply unit are included with the scope of delivery.

The connected temperature chamber(s) are addressed via this COM interface and the address that can be individually adjusted at each chamber controller.

Figure 1: Connected interface converter

4. Scope of delivery of RS485/RS422 interface converter kit

230V version

Complete kit:

Description	Art. no.
Set Interface Converter RS485/RS422 – 230V	8012-0589

Components:

Description	Art. no.
RS485/RS422 interface converter with original description of producer W & T (only needed in case of problems during start-up), including 230V power supply unit	5021-0020
2 connection cables 5m for connecting the interface converter to an individual chamber or to a plug distributor RS422 from BINDER	5023-0111
Manual for connecting the RS485/RS422 interface converter	7001-0332

115V version

Complete kit:

Description	Art. no.
Set Interface Converter RS485/RS422 – 115V	8012-0599

Components:

Description	Art. no.
RS485/RS422 interface converter with original description of producer W & T (only needed in case of problems during start-up) including 115V power supply unit	5021-0031
2 connection cables 5m for connecting the interface converter to an individual chamber or to a plug distributor RS422 from BINDER	5023-0111
Manual for connecting the RS485/RS422 interface converter	7001-0332

5. Networking

RS485 temperature chambers are connected via the RS422/RS485 interface converter.

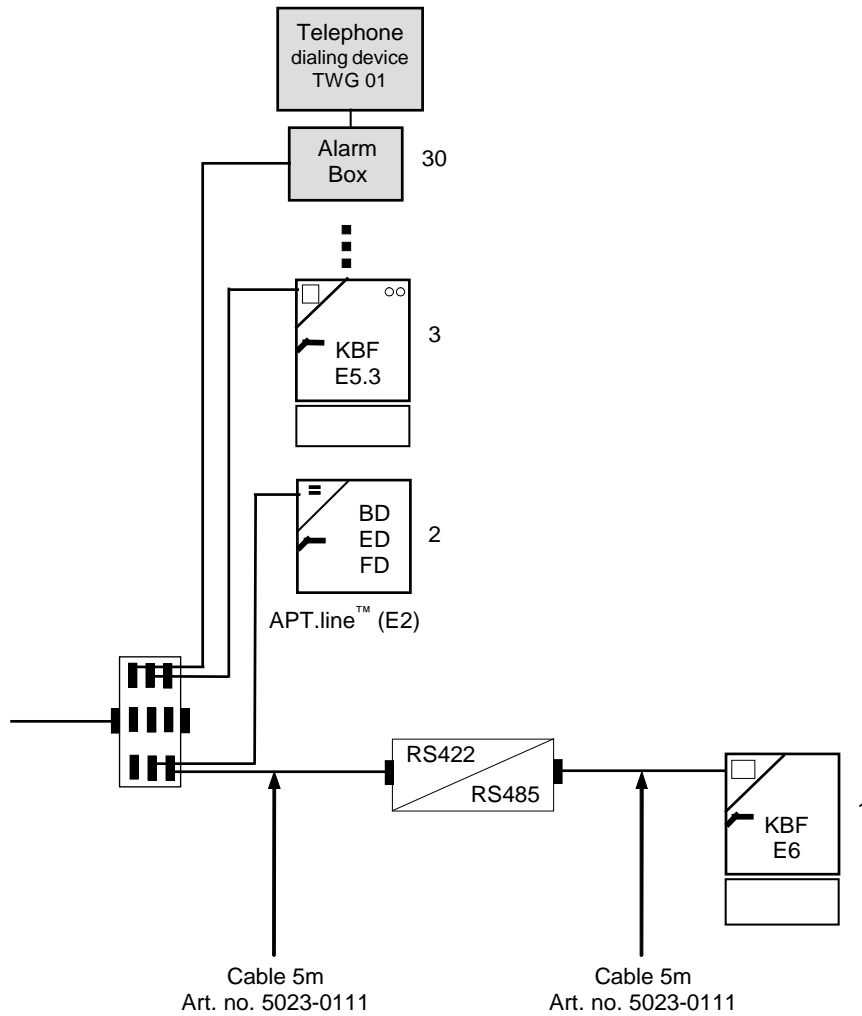
Figure 2: Networking via RS422/RS485 interface converter

The chamber address required for the unambiguous chamber identification must be entered by hand individually and only once at the controller of each temperature chamber or measurement device. For a description see the respective operating manual of the chamber or device bearing an RS485 interface.

5.1. Bus termination

To connect temperature chambers from BINDER, the recommended operating mode of the interface converter is termination enabled (BINDER factory setting).

Some controllers, like RD3, MB1, MB2, have an internal termination. When combining different controllers, communication errors can occur which are indicated by APT-COM™ as „timeout“ or „CRC error“. In this case it can be necessary to disable termination at the interface converter.

To disable termination, set DIP switches 6 and 7 in the interface converter to Off. You can safely try the correct setting. No damage of controller or interface converter will occur.

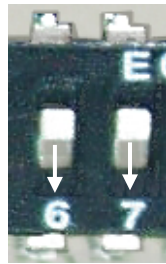
Figure 3: Termination setting via DIP switches 6 and 7

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