**Safety Instructions / 安全提示 / Инструкции за безопасност**  
**Bezpečnostní pokyny / Sicherheitshinweise / Sikkerhedshenvisninger**  
**Ohutusjuhised / Indicaciones de seguridad / Consignes de sécurité**  
**Υποδείξεις για την ασφάλεια / Sigurnosna napomena**  
**Biztonsági utasítások / Treoracha sábháilteachta / Avvertenze di sicurezza**  
**Saugumo nuorodos / Drošības norādījumi / Noti dwar is-sigurtà**  
**Veiligheidswaarschuwingen / Wskazówki bezpieczeństwa**  
**Instruções de segurança / Reguli de siguranță / Säkerhetsanvisningar**  
**Turvallisuushjeet / Bezpečnostné pokyny / Varnostna opozorila**

**KB (E7) • KB PRO (E7) • KBF (E7) • KBF PRO (E7)**

[www.binder-world.com/manuals](http://www.binder-world.com/manuals)

**Download the full operating manual.**

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---

## **KB / KB PRO (E7)**

Cooling Incubators with compressor technology and program control

## **KBF / KBF PRO (E7)**

Constant climate chambers with program control

---

### **Important note**

The present document is only valid in conjunction with the operating manual that correspond to your device

- |                                                                  |                    |
|------------------------------------------------------------------|--------------------|
| • KB (E7)                                                        | art. no. 7001-0567 |
| • KB PRO (E7)                                                    | art. no. 7001-0568 |
| • KBF + KBF PRO (E7)                                             | art. no. 7001-0569 |
| • Additionally with the ICH Q1B Light Module accessory           | art. no. 7001-0577 |
| • Additionally with the CO <sub>2</sub> Control Module accessory | art. no. 7001-0576 |

The initial commissioning and usage of the aforementioned devices (during the entire product life cycle) is only permitted with the operating manual. The operating manual must be read completely before usage.

The present document is not a replacement for the operating manual, as it only contains selected information.

**You can download the appropriate operating instructions for your device using this QR code or at [www.binder-world.com/manuals](http://www.binder-world.com/manuals). If this is not possible, please contact BINDER GmbH:**

|                 |                                                                |
|-----------------|----------------------------------------------------------------|
| Address         | Post office box 102, 78502 Tuttlingen, Germany                 |
| Tel.            | +49 7462 2005 0                                                |
| Internet        | <a href="http://www.binder-world.com">www.binder-world.com</a> |
| Service Hotline | +49 7462 2005 555                                              |
| Service Fax     | +49 7462 2005 93 555                                           |

## Safety instructions

The following represents the hazards for people. Safety information with lower risk levels, and important detailed information about safety, can be found in the operating manual. To avoid injuries and damage observe all safety instructions of the operating manual. It is very important that the operating manual is read completely before initial commissioning and usage.

### Pictograms

|                                                                                                                                                              |                                                                                                                        |                                                                                                                     |                                                                                                                                         |
|--------------------------------------------------------------------------------------------------------------------------------------------------------------|------------------------------------------------------------------------------------------------------------------------|---------------------------------------------------------------------------------------------------------------------|-----------------------------------------------------------------------------------------------------------------------------------------|
| <br>Risk of injury. Observe the safety instructions in the operating manual | <br>Electrical hazard                 | <br>Hot surface                   | <br>Explosive atmosphere                             |
| <br>Stability hazard                                                        | <br>Lifting hazard                    | <br>Scalding hazard               | <br>Risk of chemical burns                           |
| <br>Harmful substances                                                      | <br>Biohazard                         | <br>UV light hazard               | <br>CO <sub>2</sub> suffocation and poisoning hazard |
| <br>Gas cylinders                                                         | <br>Mandatory regulation            | <br>Read operating instructions | <br>Disconnect the power plug                      |
| <br>Lift with several persons                                             | <br>Lift with mechanical assistance | <br>Wear protective gloves      | <br>Wear safety goggles                            |
| <br>Do NOT touch                                                          | <br>Do NOT spray with water         | <br>Do NOT climb                |                                                                                                                                         |

|                                                                                     |                                                                                                                                                                                                                                                                                                                                                                                                                 |
|-------------------------------------------------------------------------------------|-----------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------|
|  |  <b>DANGER</b>                                                                                                                                                                                                                                                                                                               |
|  | <p><b>Electrical hazard during live maintenance work.</b><br/> <b>Deadly electric shock.</b></p> <p>Ø Do NOT remove the rear panel of the chamber.</p> <p>➤ Disconnect the chamber before conducting maintenance work. Turn off the main power switch and pull the power plug.</p> <p>➤ Make sure that general maintenance work will be conducted by licensed electricians or experts authorized by BINDER.</p> |

|                                                                                     |                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                             |
|-------------------------------------------------------------------------------------|-------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------|
|    | <div data-bbox="802 241 1061 309">  <b>DANGER</b> </div> <p><b>Explosion or implosion hazard through the introduction of unsuitable loading material, or through combustible dusts or explosive mixtures in the vicinity of the chamber or in the chamber.</b></p> <p><b>Serious injury or death from burns and / or explosion pressure.</b></p> <ul style="list-style-type: none"> <li>Ø Do NOT operate the chamber in potentially explosive areas.</li> <li>Ø Make sure that there are NO combustible dust or air-solvent mixture in the vicinity of the chamber.</li> <li>Ø Do NOT introduce any substance combustible or explosive at working temperature into the chamber, in particular no energy sources such as batteries or lithium-ion batteries</li> <li>Ø Make sure that there are NO combustible dust or air-solvent mixture in the inner chamber.</li> </ul> |
|   | <div data-bbox="802 770 1061 837">  <b>DANGER</b> </div> <p><b>Electrical hazard by water entering the chamber.</b></p> <p><b>Deadly electric shock.</b></p> <ul style="list-style-type: none"> <li>Ø Do NOT install the chamber in damp areas or in puddles.</li> <li>➢ Set up the chamber in a splash-proof manner.</li> <li>➢ Before cleaning, turn off the chamber at the On/Off switch and disconnect the power plug. Let the chamber cool down to ambient temperature.</li> <li>Ø The chamber must NOT become wet during operation, cleaning, or maintenance. Do NOT spill water or cleaning agents over the inner and outer chamber surfaces.</li> <li>➢ Completely dry the chamber before turning it on again.</li> </ul>                                                                                                                                          |
|  | <div data-bbox="802 1267 1061 1335">  <b>DANGER</b> </div> <p><b>Electrical hazard through overload of contacts on the zero-voltage relay alarm output.</b></p> <p><b>Deadly electric shock. Damage to the switching contacts and connection socket.</b></p> <ul style="list-style-type: none"> <li>Ø Do NOT exceed the maximum switching load of 24V AC/DC – 2.5A.</li> <li>Ø Do NOT connect any devices with a higher loading capacity.</li> </ul>                                                                                                                                                                                                                                                                                                                                                                                                                     |
|  | <div data-bbox="802 1572 1061 1639">  <b>DANGER</b> </div> <p><b>Electrical hazard due to missing protective conductor connection or high leakage current in case of improper installation.</b></p> <p><b>Deadly electric shock.</b></p> <ul style="list-style-type: none"> <li>➢ Make sure that the chamber's power plug and the power socket match and securely connect the electrical protective conductors of the chamber and the house installation.</li> <li>➢ Only connect the plug to a socket if it has an intact earth connection.</li> </ul>                                                                                                                                                                                                                                                                                                                  |

|                                                                                                                                                                     |                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                   |
|---------------------------------------------------------------------------------------------------------------------------------------------------------------------|---------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------|
|   | <div style="background-color: #ffcc00; text-align: center; padding: 5px;">  <b>WARNING</b> </div> <p><b>Danger of intoxication and infection through contamination of the chamber with toxic, infectious or radioactive substances.</b></p> <p><b>Poisoning. Damages to health.</b></p> <ul style="list-style-type: none"> <li>Ø Do NOT introduce any substance which could lead to release of toxic gases</li> <li>➤ Protect the interior of the chamber from contamination by toxic, infectious or radioactive substances.</li> <li>➤ Take suitable protective measures when introducing and removing toxic, infectious or radioactive material.</li> <li>Ø NEVER take a chamber contaminated with toxic substances or sources of infection for recycling according to Directive 2012/19/EU.</li> <li>➤ Prior to disposal, remove all toxic substances and sources of infection from the chamber.</li> <li>➤ A chamber from which all toxic substances or sources of infection cannot be safely removed must be considered as “special” waste according to national law. Dispose of it accordingly.</li> </ul> |
|---------------------------------------------------------------------------------------------------------------------------------------------------------------------|---------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------|

|                                                                                                                                                                                                                                                                                                                                               |                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                    |
|-----------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------|------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------|
|     | <div style="background-color: #ffff00; text-align: center; padding: 5px;">  <b>CAUTION</b> </div> <p><b>Risk of injury and damages by lifting heavy loads and by sliding or tilting of the chamber due to improper lifting or transport.</b></p> <p><b>Injuries.</b></p> <ul style="list-style-type: none"> <li>➤ For transport, secure the chamber with transport straps.</li> <li>➤ Lift chambers sizes 65 and 130 with the aid of 4 people, chambers size 260 with the aid of 6 people or use a fork lifter.</li> <li>➤ Lift chambers sizes 470, 720, 1060 and 1600 using technical devices (fork lifter).</li> <li>Ø Do NOT load the lower housing cover with heavy objects while the chamber door is open and do NOT climb on it.</li> </ul> |
|-----------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------|------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------|

|                                                                                     |                                                                                                                                                                                                                                                                                                                                                                                                      |
|-------------------------------------------------------------------------------------|------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------|
|  | <div style="background-color: #ffff00; text-align: center; padding: 5px;">  <b>CAUTION</b> </div> <p><b>Danger of burning by touching hot chamber parts during or after operation.</b></p> <p><b>Burns.</b></p> <p>Do NOT touch the inner surfaces, the glass doors or the loading material during operation.</p> |
|-------------------------------------------------------------------------------------|------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------|

|                                                                                                                                                                                                                                                                                                                                                 |                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                           |
|-------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------|-------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------|
|     | <div style="background-color: #ffff00; text-align: center; padding: 5px;">  <b>CAUTION</b> </div> <p><b>Danger of chemical burns through contact with skin or ingestion of the neutral cleaning agent or through eye contact with the disinfectant spray.</b></p> <p><b>Skin and eye damage.</b></p> <ul style="list-style-type: none"> <li>➤ Do not ingest the neutral cleaning agent. Keep it away from food and beverages.</li> <li>➤ Wear protective gloves and goggles.</li> <li>➤ Avoid skin contact with the neutral cleaning agent.</li> </ul> |
|-------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------|-------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------|

For chambers with the ICH Q1B Light Module accessory:

|  |  <b>WARNING</b>                                                                                                                                                                                                                                                                    |
|-----------------------------------------------------------------------------------|---------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------|
|                                                                                   | <p><b>Danger of eye and skin injury by UV light hazard.</b><br/> <b>Eye and skin injury.</b></p> <ul style="list-style-type: none"> <li>Ø Do NOT look directly into the light.</li> <li>➤ Minimize skin exposure.</li> <li>➤ Wear UVA protective goggles when opening the door of the cooling incubator / constant climate chamber when the light is on.</li> </ul> |

For chambers with the CO<sub>2</sub> Control Module accessory:

|  |  <b>DANGER</b>                                                                                                                                                                                                                                                                                                                                                             |
|-----------------------------------------------------------------------------------|-------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------|
|                                                                                   | <p><b>Danger of suffocation and poisoning by high concentration of CO<sub>2</sub> (&gt; 4 Vol.-%).</b><br/> <b>Death by suffocation.</b></p> <ul style="list-style-type: none"> <li>Ø Do NOT set up chambers in non-ventilated recesses.</li> <li>➤ Ensure technical ventilation measures.</li> <li>➤ Observe the relevant regulations for handling CO<sub>2</sub>.</li> <li>➤ Close the CO<sub>2</sub> supply when decommissioning the chamber.</li> </ul> |

|  |  <b>WARNING</b>                                                                                                                                                                                                       |
|-------------------------------------------------------------------------------------|----------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------|
|                                                                                     | <p><b>Risk of injury through sudden release of the stored pressure energy when the valve safety is torn off.</b><br/> <b>Injuries.</b></p> <ul style="list-style-type: none"> <li>➤ Secure gas cylinders against falling (chaining).</li> <li>➤ Transport gas cylinders with a cylinder cart.</li> </ul> |

|  |  <b>WARNING</b>                                                                                                                                                                                                |
|-------------------------------------------------------------------------------------|---------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------|
|                                                                                     | <p><b>Risk of injury through sudden release of the stored pressure energy when opening the cylinder valve of a not connected cylinder.</b><br/> <b>Injuries.</b></p> <ul style="list-style-type: none"> <li>➤ Close the gas cylinder valve before connecting or removing the gas hose.</li> </ul> |

---

## Intended use

---

The chambers are suitable for exact conditioning of harmless materials.

Any solvent must not be explosive and flammable. A mixture of any component of the charging material with air must NOT be explosive. The operating temperature must lie below the flash point or below the sublimation point of the charging material. Any component of the charging material must NOT be able to release toxic gases. Other applications are NOT approved.

The chambers are designed for setting up inside a building (indoor use). The requirements described in the Operating Manual for installation site and ambient conditions must be met

The chambers are not classified as medical devices as defined by Regulation (EU) No 2017/745.

The loading material shall not contain any corrosive ingredients that may damage the machine components made of stainless steel, aluminum, and copper. Such ingredients include in particular acids and halides. Contamination of the chamber by toxic, infectious or radioactive substances must be prevented. The chambers do not dispose of any measures of explosion protection.

Details can be found in the operating manual in the chapter "Intended use". Any use of the chambers that does not comply with the requirements specified in the Operating Manual shall be considered improper use. Make sure that you have the Operating Manual that match your chamber, including any accessories.

You may not make any modifications to the chamber or device yourself, as these may limit proper usage.

To operate the chamber, use only original BINDER accessories or accessories / components from third-party suppliers authorized by BINDER. The user is responsible for any risk arising from using unauthorized accessories

BINDER GmbH is responsible for the safety features of the chamber only, provided skilled electricians or qualified personnel authorized by BINDER perform all maintenance and repair, and if components relating to chamber safety are replaced in the event of failure with original spare parts. The user is responsible for any risks arising from using unauthorized accessories/components.

### For KBF/KBF PRO:

#### Types of suitable water quality for freshwater supply:

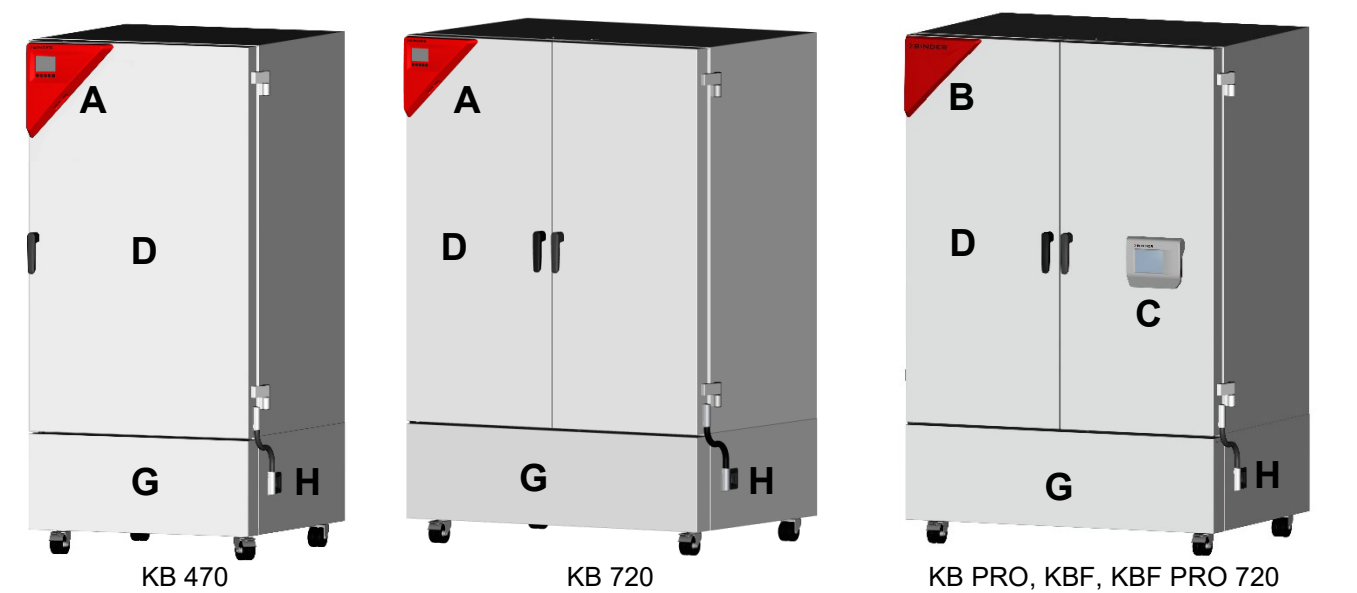
- Deionized water from a water treatment installation already existing at the customer's site. Conductivity from 1  $\mu\text{S}/\text{cm}$  up to a maximum of 20  $\mu\text{S}/\text{cm}$ . (Water, which is in equilibrium with the  $\text{CO}_2$  in the air, and has a conductivity below 1  $\mu\text{S}/\text{cm}$  (ultrapure water), may cause acid corrosion due to its low pH)
- Water treated by the optional water treatment system BINDER Pure Aqua Service (disposable system).

### For chambers with the $\text{CO}_2$ Control Module accessory:

The chambers allow the setting of an exact  $\text{CO}_2$  atmosphere inside. The chambers are not completely gas-tight. Vent out any  $\text{CO}_2$  gas that may escape via good room ventilation or a suitable connection to an exhaust system. The installation of a  $\text{CO}_2$  warning system is mandatory.

Chamber overview

Front

- A

Triangular instrument panel  
(KB all sizes, KB PRO / KBF/ KBF PRO sizes 130, 260)
- B

Triangular signal panel (KB PRO / KBF/ KBF PRO starting with size 470)
- C

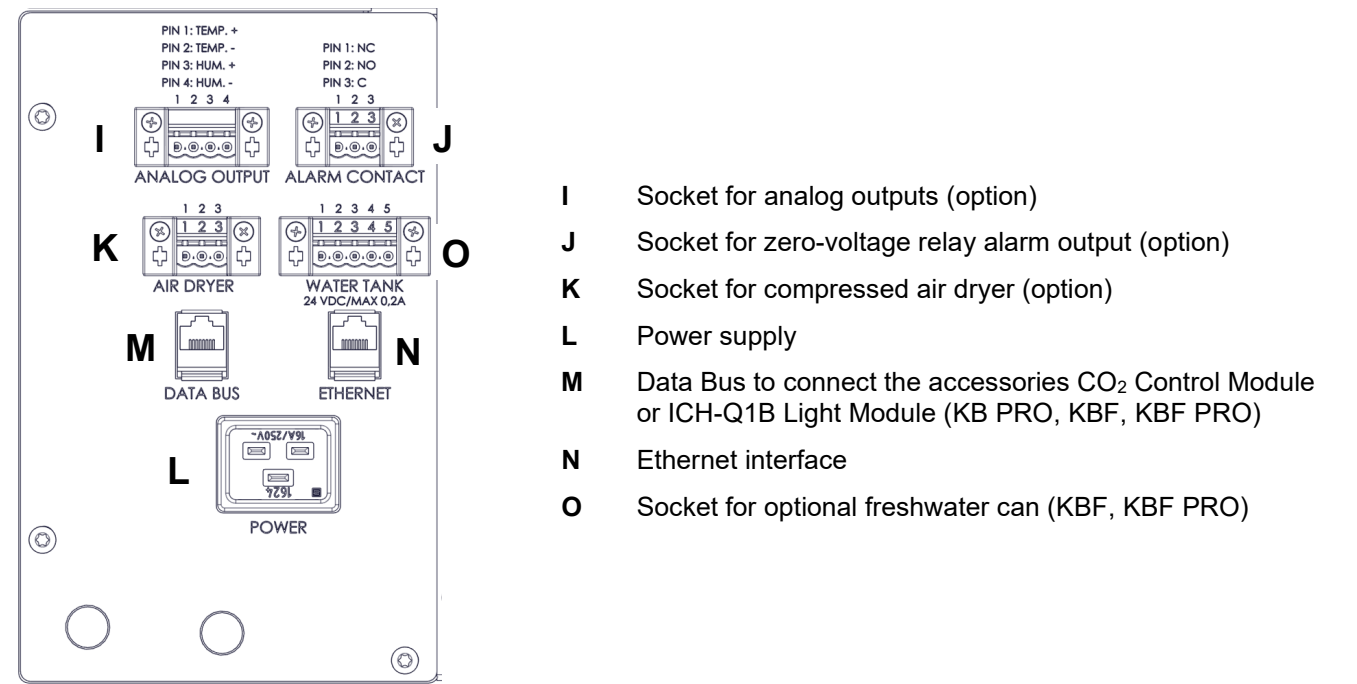
Instrument panel (KB PRO / KBF/ KBF PRO starting with size 470)
- E

Outer door(s)
- G

Technical installation room cover
- H

On/Off switch

Connection panel on the chamber rear

## Electrical connection data and installation

- ▶ The chambers come with an IEC connector plug with a power connection cable of 2 m in length (optional 5 m) and a grounded plug. Only use original connection cables from BINDER according to the specification given in the Operating Manual.
- ▶ Pollution degree acc. to IEC 61010-1: 2. Over-voltage category acc. to IEC 61010-1: II. Housing protection acc. to EN 60529: IP 20
- ▶ KB, KBF: 220-240 V Nominal voltage (+/-10%) at 50 and 60 Hz, Current type 1N~  
KB PRO, KBF PRO: 208-240 V Nominal voltage (+/-10%) at 50 and 60 Hz, Current type 1N~  
KB-UL, KBF-UL: 120 V Nominal voltage (+/-10%) at 50 and 60 Hz, Current type 1N~
- ▶ Internal over-current release category B 2 poles 16 Amp
- ▶ Nominal power

| Model, Size |    | 65  | 130  | 260  | 470  | 720  | 1060 | 1600 |
|-------------|----|-----|------|------|------|------|------|------|
| KB          | kW | 0.9 | 1.2  | 1.3  | 1.3  | 1.5  | 1.5  | ---  |
|             | A  | 3,8 | 5.0  | 5.4  | 5.4  | 6.3  | 6.3  | ---  |
| KB-UL       | kW | 0.9 | 1,2  | 1.3  | 1.3  | 1.5  | 1,5  | ---  |
|             | A  | 7.5 | 10.0 | 10.8 | 10.8 | 12.6 | 12.6 | ---  |
| KB PRO      | kW | --- | 1.0  | 1.5  | 1.5  | 1.7  | 1.8  | 1.8  |
|             | A  | --- | 4.2  | 6.3  | 6.3  | 7.1  | 7.5  | 7.5  |
| KBF         | kW | --- | 1.6  | 1.4  | 1.4  | 1.6  | 1.6  | 1.7  |
|             | A  | --- | 6.7  | 5.8  | 5.8  | 6.7  | 6.7  | 7.1  |
| KBF-UL      | kW | --- | 1.6  | 1.4  | 1.4  | 1.6  | 1.6  | 1.7  |
|             | A  | --- | 13.4 | 11.6 | 11.6 | 13.4 | 13.4 | 14.2 |
| KBF PRO     | kW | --- | 1.9  | 1.6  | 1.6  | 1.8  | 1.9  | 1.9  |
|             | A  | --- | 7.9  | 6.7  | 6.7  | 7.5  | 7.9  | 7.9  |

- ▶ The domestic socket must also provide a protective conductor. Make sure that the connection of the protective conductor of the domestic installations to the chamber's protective conductor meets the latest technology. The protective conductors of the socket and plug must be compatible!
- ▶ When connecting, please observe the regulations specified by the local electricity supply company as well as the local or national electrical regulations.
- ▶ Observe a sufficient current protection according to the number of devices that you want to operate. We recommend the use of a residual current circuit breaker.
- ▶ Prior to connection and start-up, check the power supply voltage. Compare the values to the specified data located on the chamber's type plate.
- ▶ To completely separate the chamber from the power supply, you must disconnect the power plug. Install the chamber in a way that the power plug is easily accessible and can be easily pulled in case of danger.

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